

Celloger[®] Nano

Benchtop digital microscope

| Quick Manual




red dot winner 2022

Quick Installation

1. Hardware

1.1 Wireless connection

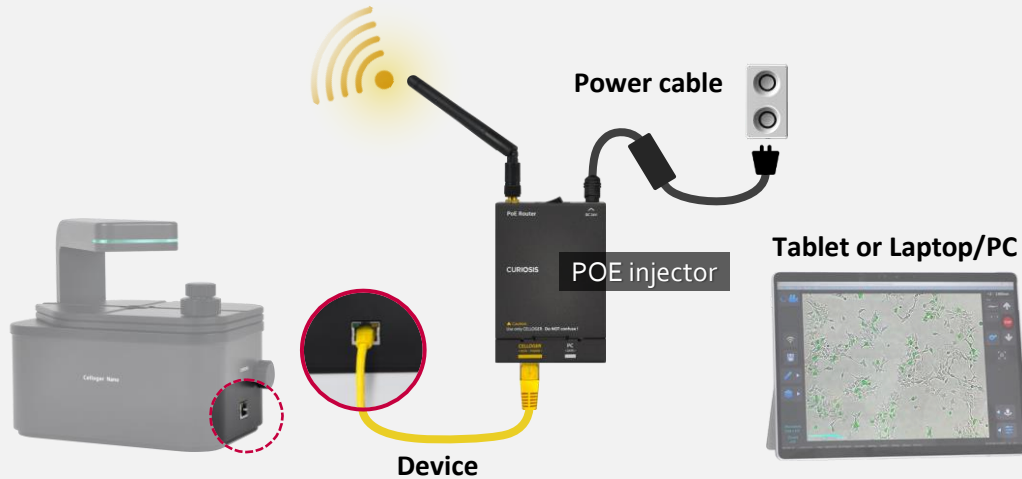


Step 1. Connect the **LAN cable** and **Power cable** to the **POE injector**.

Step 2. Connect the **LAN cable** to the **device**.

Step 3. Connect the **Power cable** to an **outlet**.

1.2 Wired connection

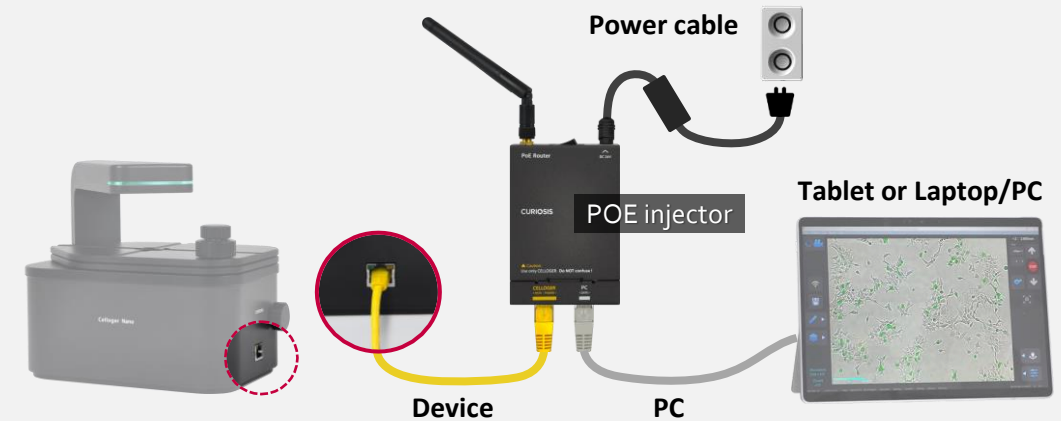

Step 1. Connect **two LAN cables** and a **Power cable** to the **POE injector**.

Step 2. Connect the **LAN cable(yellow)** to the **device** and another **LAN cable(white)** to the **PC**.


Step 3. Connect the **Power cable** to an **outlet**.




The device can be connected to **Windows-based(more than Window 10 and above)** tablets, laptops, and PCs.

Quick Installation


2. Software

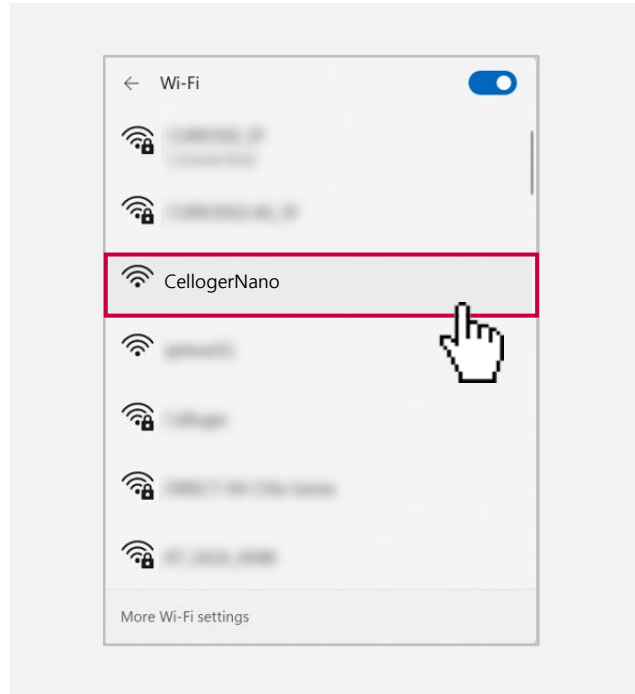

Step 1. Go to the Wi-Fi settings of the tablet and connect to the network named “**CellogerNano**”.

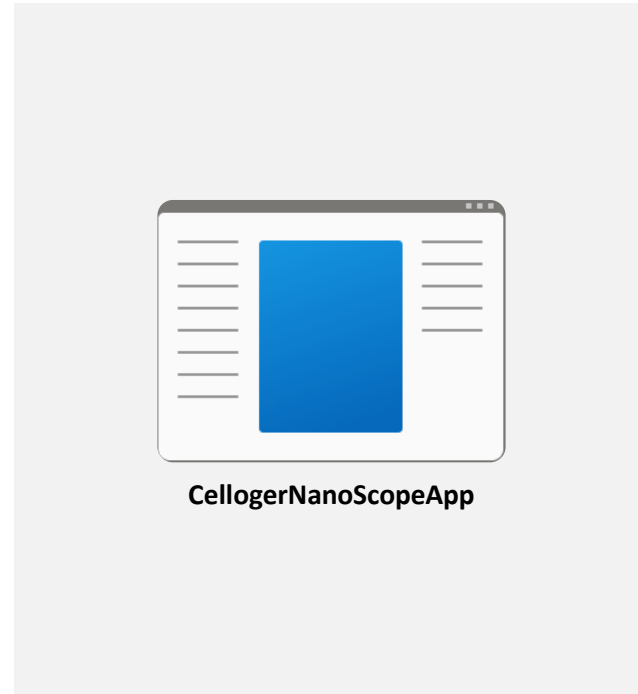

Step 2. Open **CellogerNanoScopeApp**.

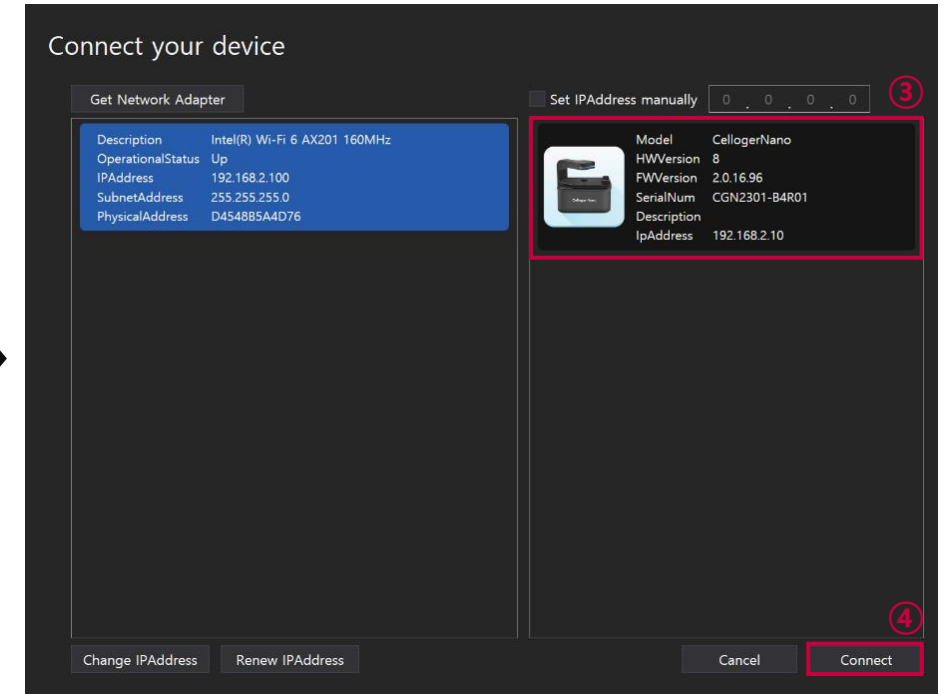

Step 3. Press the **device icon**.

Step 4. Press the **Connect** to connect the device.

***Note:** If the device icon does not appear when using a wired connection, please refer to the **full manual** for instructions on adjusting the PC network settings.

Thank you

End of Document

Curiosis Inc.

4F, 10, Teheran-ro 38-gil, Gangnam-gu, Seoul 06221, South Korea | T 82 2 508 5237 | F 82 2 508 5246 | sales@curiosis.com
www.curiosis.com